# DESIGNBOOK

## LABORATORY INNOVATION

**BICASA** 

 BICASA SRL  
Bernareggio  
Italy

website  
[WWW.BICASA.IT](http://WWW.BICASA.IT)

**BICASA** 

---

**GROUP**

---

## SOMMARIO / INDEX

<b>LA NOSTRA AZIENDA</b> Our company	4
<b>IDEE PER LABORATORIO</b> Laboratory concept	8
<b>GESTIONE DEL PROGETTO</b> Project management	12
<b>BANCHI DA LABORATORIO MECC</b> Mecc laboratory benches	16
<b>PARETE TECNICA DOCKWALL</b> Dockwall partition system	26
<b>CAPPE CHIMICHE RX</b> RX fume cupboards	32
<b>IKARO &amp; KOMPATTO</b> Supply beams & mobile workstations	48
<b>MODULI VERTICALI COLUMNIA</b> Columnia service column	64
<b>KRISTAL VSM</b> Vertical service module	70
<b>ARMADI DI SICUREZZA</b> Storage safety cabinets	74
<b>CONTENITORI PER LABORATORIO</b> Laboratory cabinets	78
<b>PIANI DI LAVORO</b> Worktops	88

# LA NOSTRA AZIENDA

OUR COMPANY

## LA STORIA / HISTORY

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*Settled in Sesto S. Giovanni (Milan), Bicasa starts its own rise in 1947 from the initiative of the two founders, Luigi Biffi and Giuseppe Casiraghi. Postwar economic boom allows the company to suddenly become a leading supplier for laboratory market. Following a constant innovation, Bicasa becomes a successful enterprise in glassware processing (blowing), **scientific instruments manufacturing and technical furniture designing**, employing up to 100 human resources with 18.000 produced items.*

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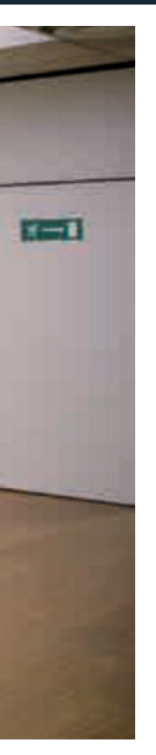
*Tradition and experience have been knowing a permanent renewal till third generation, making BICASA a leading supplier in italian market for laboratory furniture. The international markets presence is assured by an open minded view of globalization, focusing on foreign markets needs.*

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*Maintaining and expanding our position as a leading supplier in the worldwide laboratory furniture market. We are focused on selection and training of highly specialized and motivated employees and distribution partners dedicated to provide professional and excellent consulting to our customers. We are strongly committed to achieve our goals based on the **guiding lines of honesty, trust, respect and quality**.*

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*The headquarter office is in Bernareggio, close to Milan. The building, placed on 20.000 sqm overall, includes offices and manufacturing departments for wood and steel processing. Design and production are wholly factory-made through dedicated internal divisions, ensuring a total flexibility of manufacturing cycle.*

## SISTEMA DI QUALITA' CERTIFICATO ISO 9001 ISO 9001 QUALITY SYSTEM CERTIFICATION

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*Since 1996 BICASA **works in accordance with ISO 9001 quality management system**, for quality assurance in design, development, production, installation and servicing.*

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*Equipped benches, partition systems, supply beams, fume cupboards and furnishing complements set up the **integrated system** of BICASA collection, designed to offer a **complete solution** for all laboratory needs.*

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*Constant development of instruments and new methodologies applied in scientific research, set up Bicasa guidelines to **get technological innovations** for multipurpose furnishing systems.*

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*Raw materials selection, chromatic research and attention to details are expressed through **stackable elements**, following a furniture concept which design and functionality join to get a unique output in.*

## QUALITA' E SICUREZZA / QUALITY AND SAFETY STANDARDS

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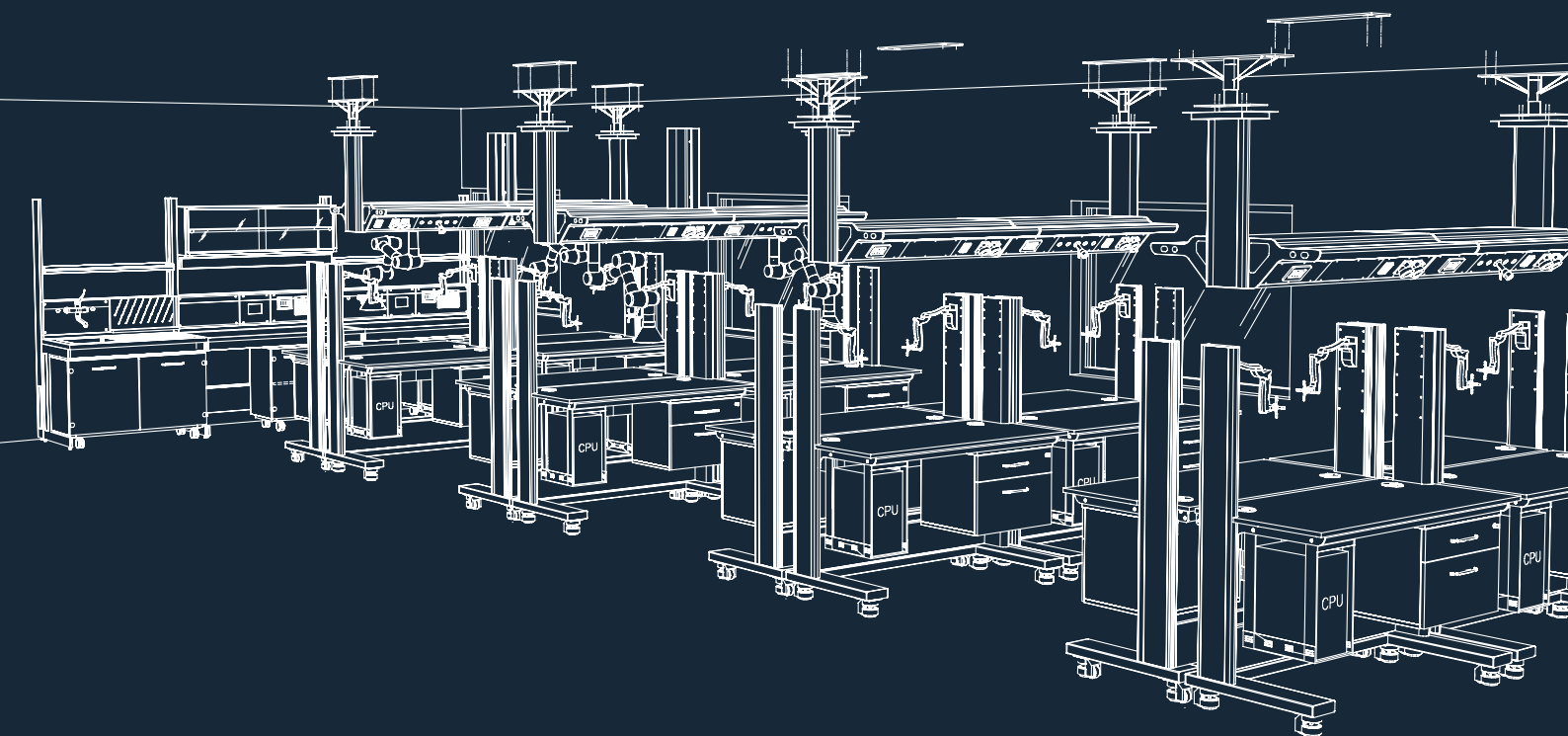
*Manufacturing quality and safety standards, fitted with research and design, are the mission of a company always focused on customers satisfaction. Furniture systems are tested and certified by third parties, assuring the **highest product quality and the strictest safety standards observance** for laboratory environments.*

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*Materials and finishes involved in furniture production have specific laboratory features to join together **chemical-mechanical resistance and surfaces decontamination requirements**. Selection of fireproof parts, matched with technological services and fittings according to international standards, are warranties for operative safety.*

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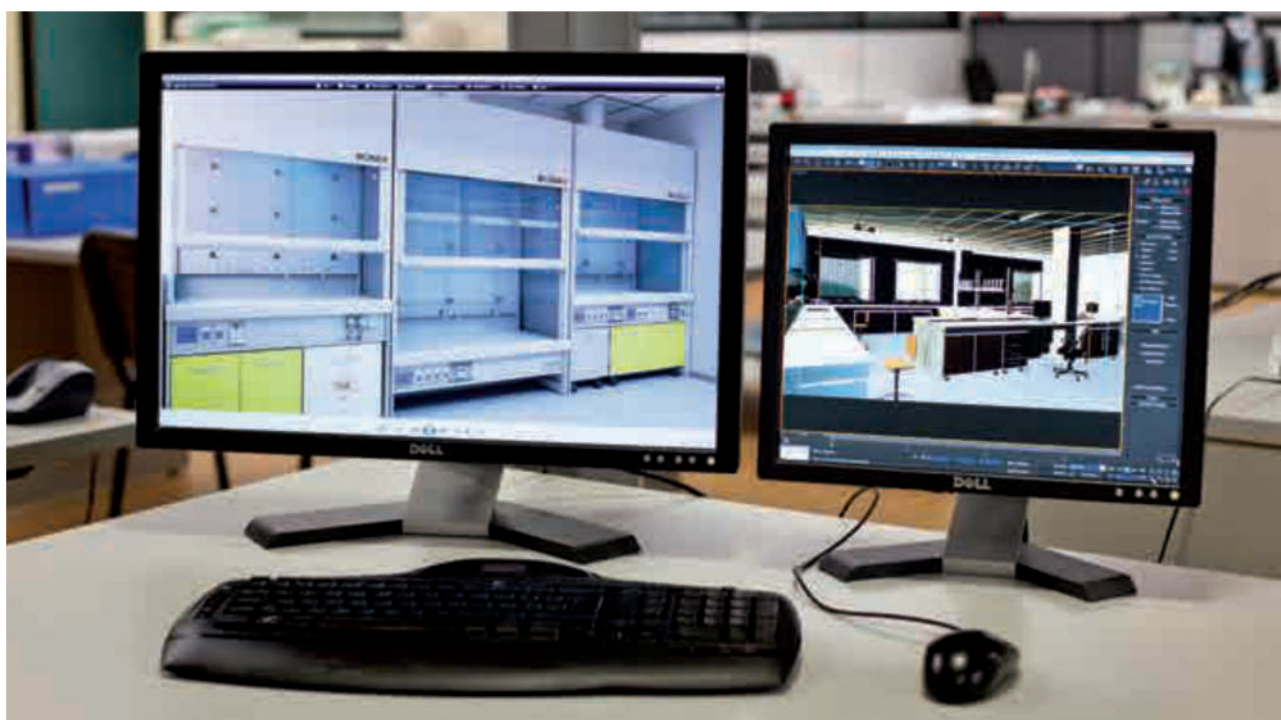
*Specific requirements evaluation, improvement advises and customized solutions for each project are the strength points that characterize Bicasa, not just as products supplier, but **problems solving partner**.*

## KNOW-HOW TECNICO / TECHNICAL KNOW-HOW

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*The experience gained since 1947 has given high skills of comprehension in laboratories processes to the company. Bicasa offers **complete and advanced know-how** to support customer operative requirements.*

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*During the project planning phase each individual element is analysed in relation to existing rooms and intended use. The **development of CAD 2D - 3D drawings and rendering presentations** make the customer able to define single details of execution.*

## ASSISTENZA IN LOCO / ON SITE MANAGEMENT

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*Really often projects require integration with existing or in process architectural works. Bicasa offers a complete **synergy work with architects and construction firms**, achieving a perfect integrated performance.*

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*Laboratory benches are certified according to **EN 13150**, European standard concerning “Working benches for laboratories: dimensions, safety requirements and test methods”.*

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*Laboratory benches are certified according to **BGR234/09.06** German standard, concerning workplace safety for storage devices. The load capacity tested and certified is up to 400 kgs uniformly distributed, for all workstation types.*

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*A modern lab requires highly adaptable systems able to provide integration with hi-tech equipment, making possible for end-users to operate in an ergonomic and functional workplace. MECC furniture system is designed to rationalize spaces, to distribute supply lines and to place service outlets or accessories on the most comfortable position. Through **the variable combination of each element** it is possible to arrange the configuration required.*

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***Aluminum profiles**, used to design service spine vertical uprights, ensure a stronger structural rigidity thanks to the higher mechanical precision of the assembly and reduce the overall weight of MECC benches. Aluminum, a recyclable material, improves significantly the resistance to corrosion, even in case of abrasion of the polyurethane powder coating.*

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*Bench service spine is designed to accept service outlets and supply lines distribution, placed in a protected and hidden compartment with an overall depth of 150 mm. Side uprights make bench able to be arranged with variable vertical combinations and install overhead cabinets or shelves, in order to best rationalize space use. Modular elements can be horizontally connected to reach the dimension required.*

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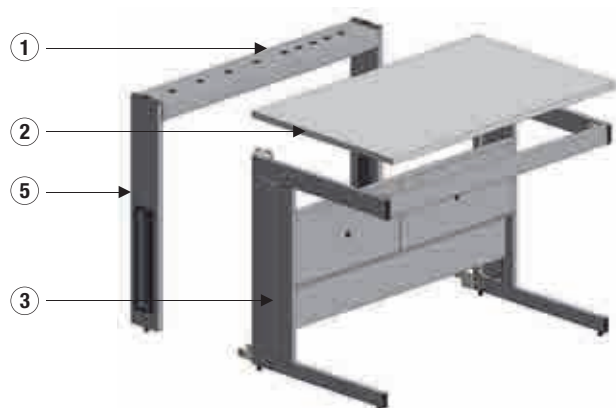
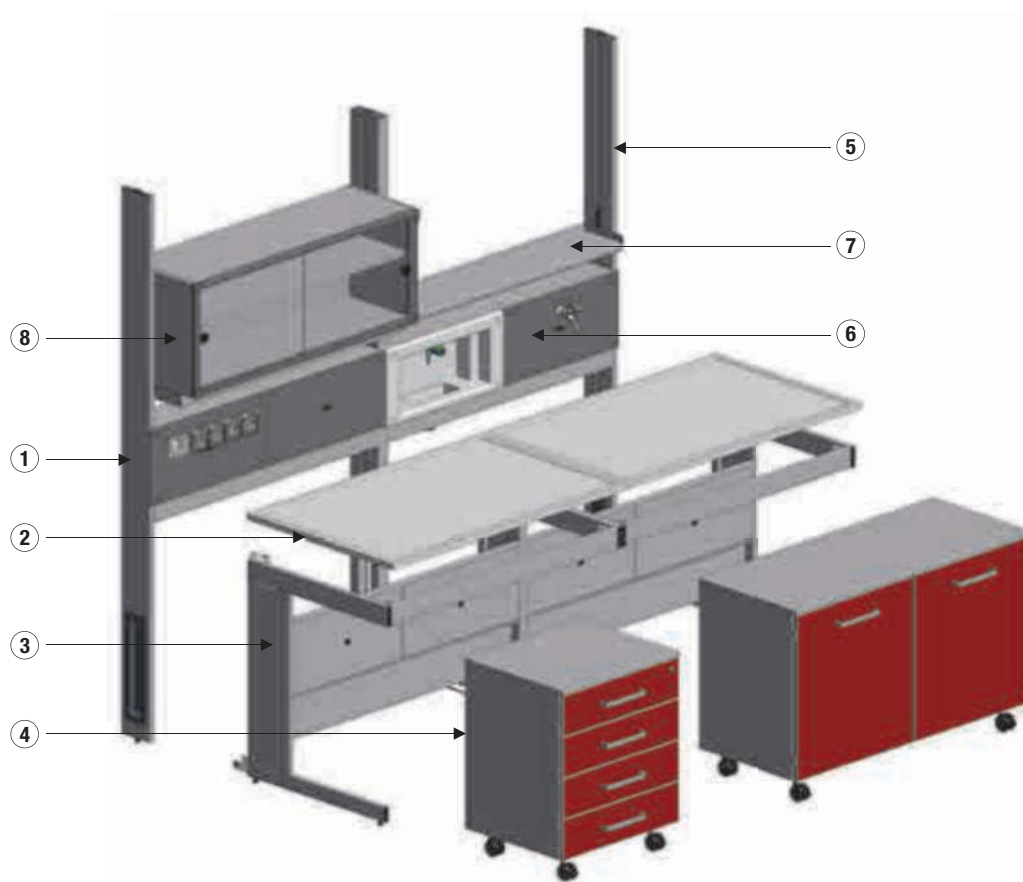
*Technical panels are designed to provide a functional distribution of service outlets, easy to be upgraded and changed. **The horizontal combination of modular panels can be also vertically extended** with additional service levels, in order to install a great number of service outlets. Panels are made of corrosion resistant molded plastic and can accept any kind of service outlet, built-in cupsink unit included.*

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*The partition system is designed in compliance with the requirements of laboratory use, applying dedicated solutions with materials suitable for the intended use. Supply lines distribution and service outlets arrangement are fully integrated in a close compartment, with an access panel - removable even if the workstation is installed. The aluminium vertical uprights, provided with a rail for suspended accessories, are ready to accept height adjustable shelves and/or overhead cabinets.*

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*Masonry walls have structural limits due to supply lines or service outlets installation. The laboratory partition system makes possible to integrate every element required, with a modular frame easy to be rearranged and fully accessible for maintenance.*

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*The partition wall consists of aluminium uprights and crossbars, ready to accept the vertical closing elements. The construction of the filler panels can be made either of powder coated steel or resin-bonded particleboard. The flexibility of the system makes possible to include frame with double glazing (laminated safety glass), hinged or sliding doors and corner junctions between perpendicular walls.*

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*The partition wall has a total thickness of mm 200, designed to accept in a dedicated compartment all the supply lines, drainage system included. The holes prearranged on the aluminium profiles provide the possibility to install the fittings with a vertical and horizontal distribution up to the final point-of-use. The plastic made modular service panels make available functional arrangements of service outlets, easy to be changed.*

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
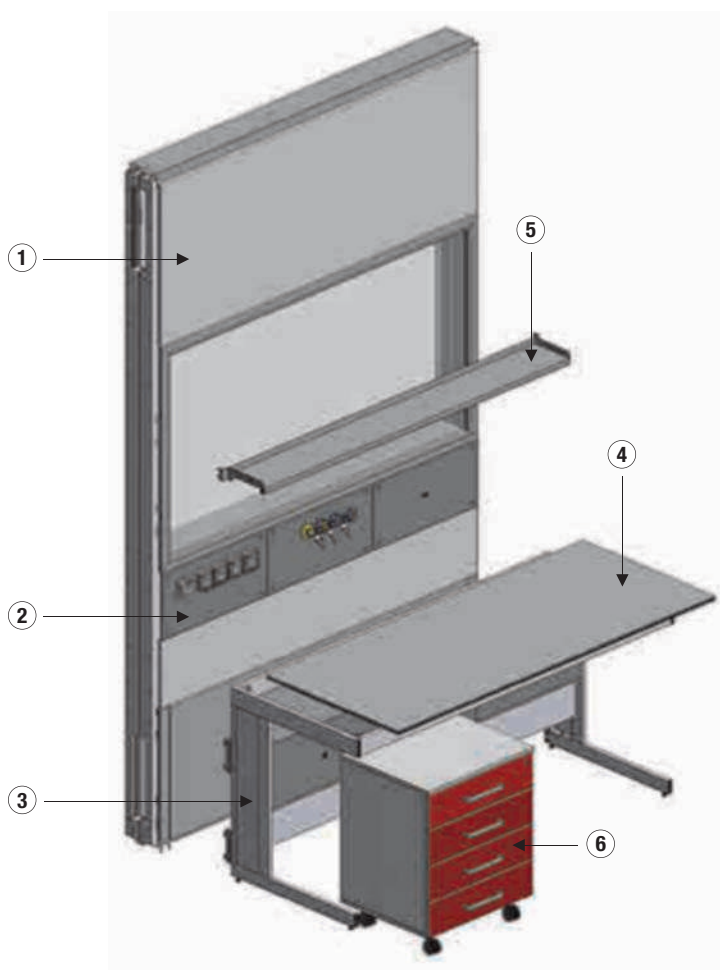
*The partition wall is a system consisting of single elements joined together in order to reach the dimension required. The modular construction, with a wide range of width measures available, provides the possibility to install laboratory tables or mobile workstation with equal dimensions to achieve a perfect combination.*

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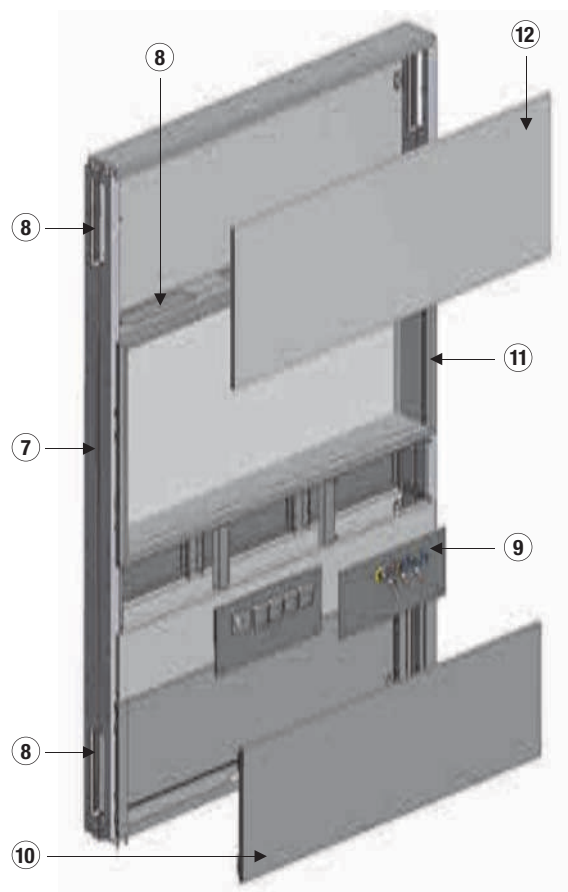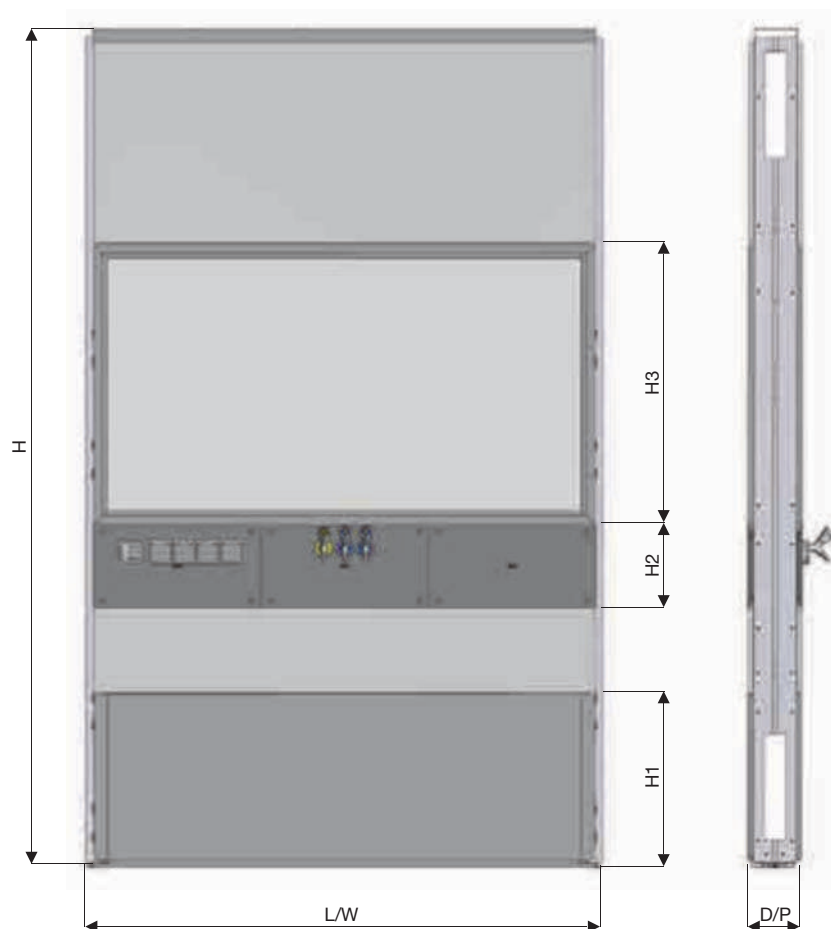
*Since 1947, Bicasa has been gaining a solid experience in laboratory design so that the different requirements of end users are fully understood by a highly skilled team of people. Safety in hazardous working conditions is a must for the design and construction of protection equipment. RX fume cupboards represent the achievement of our continuous effort as far as high-quality materials, innovative technical solutions and high performances for risk containment are concerned.*

## MATERIALI HI-TECH / HIGH-TECH MATERIALS

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*Framework is fully manufactured in aluminium - 100% recyclable material - consisting of Bicasa custom-made extruded profiles. The round-shaped design is the distinctive feature of RX fume cupboards, which improves end user comfort and increases airflow efficiency.*

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*The wide side glasses, available on RX SMART, improve the workspace visibility and make the end user able to check the processes inside the fume cupboard. The panels, made of safety laminated glass, are fixed to the frame using snap-fit aluminium profiles.*

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*Mechanical parts, operating in the vertical sliding movement, are installed in an external compartment, preventing degradation due to corrosive agents inside the working chamber. In case of breakage of either one or both belts simultaneously, the safety block system of the independent counterweights prevents sash from falling down.*

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*RX fume cupboards are a dynamic and modular system, to be easily customized in compliance with end user requirements. RX SMART, with its slim side walls, offers more internal space available and also side glasses to increase the workspace illuminance. RX FULL is designed with vertically arranged services, directly installed on side wall modules and ready to accept a great number of utilities, in order to increase flexibility of the workspace.*

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*RX is designed to offer maximum performances and safety to end users. Fume hoods are tested and certified with face velocity values of 0.15 m/s, 0.30 m/s and 0.50 m/s, to be used depending on working conditions. All tests reported excellent results of containment and robustness without any supportive airflow.*

<b>FACE VELOCITIES:</b>	0.15 - 0.30 - 0.50 m/s	testate e certificate / <i>tested and certified</i>
<b>CONTAINMENT:</b>	≤ 0.01 ppm	tutti i modelli e velocità / <i>all models and face velocities</i>
<b>ROBUSTNESS:</b>	≤ 0.298 ppm	Valore massimo con velocità frontale di 0.15 m/s per cappe con camera fino a mm 1700 / <i>Maximum value achieved with a 0.15 m/s face velocity and workspace width up to mm 1700</i>

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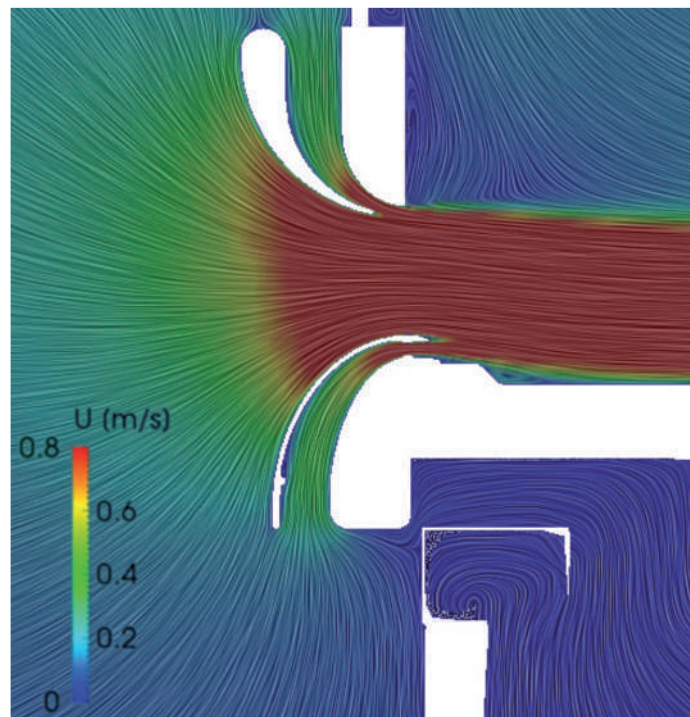
*RX fume cupboards are certified according to **EN 14175 standard Part 2-3** “Laboratory fume cupboards” - **Part 6** “Variable air volume fume cupboards” and **ASHRAE 110-2016**.*

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*The research applied to airflows entering the workspace and the construction of internal compartment are the reasons of achievement of such high performances. In order to achieve the best technical solutions, the CFD analysis (Computational Fluid Dynamics), calculated on RX mathematic model, allowed accurate evaluations under different conditions of use.*

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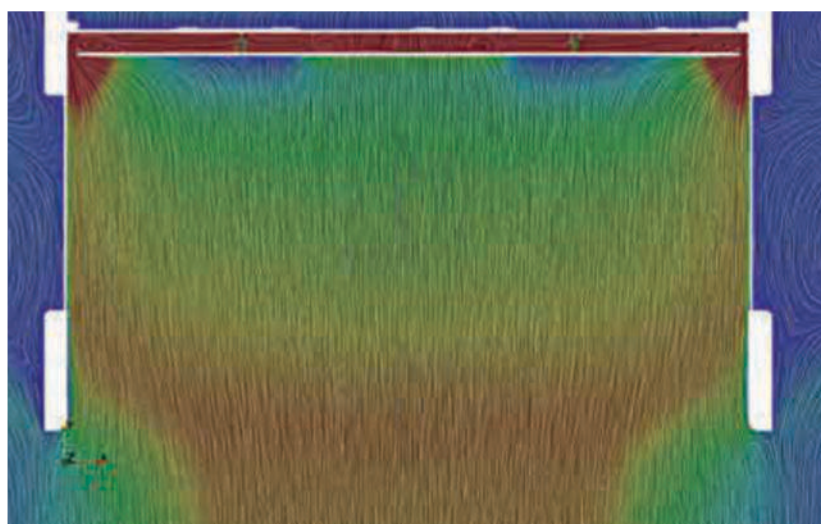
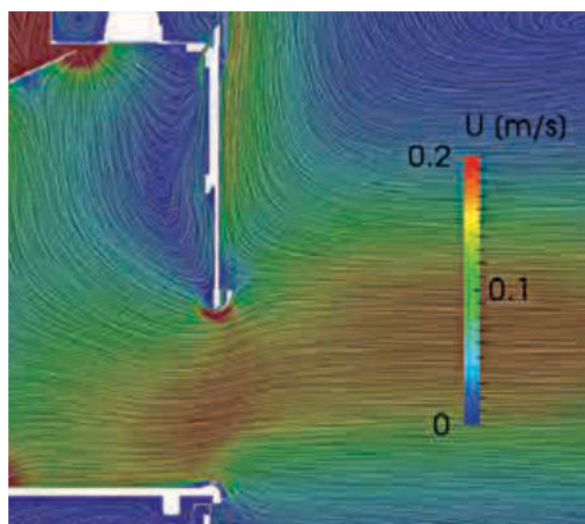
*The rated power of the equipment does not include the power consumption of sockets and remote centrifugal fan.*

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
Air volume flow rate tested and certified with face velocity equal to V1=0.15 m/s, V2=0.30 m/s, V3=0.50 m/s

CAPPE CHIMICHE WALK IN WALK IN FUME CUPBOARDS			RX FULL 1.5WI	RX FULL 1.8 WI	RX FULL 2.1WI	RX FULL 2.4WI
PORTATA ARIA ESTRATTA AIR VOLUME FLOW RATE		m³/h	1220	1490	1830	2200
CONTENIMENTO INNER-PLANE INNER-PLANE CONTAINMENT		ppm	≤ 0.01	≤ 0.01	≤ 0.01	≤ 0.01

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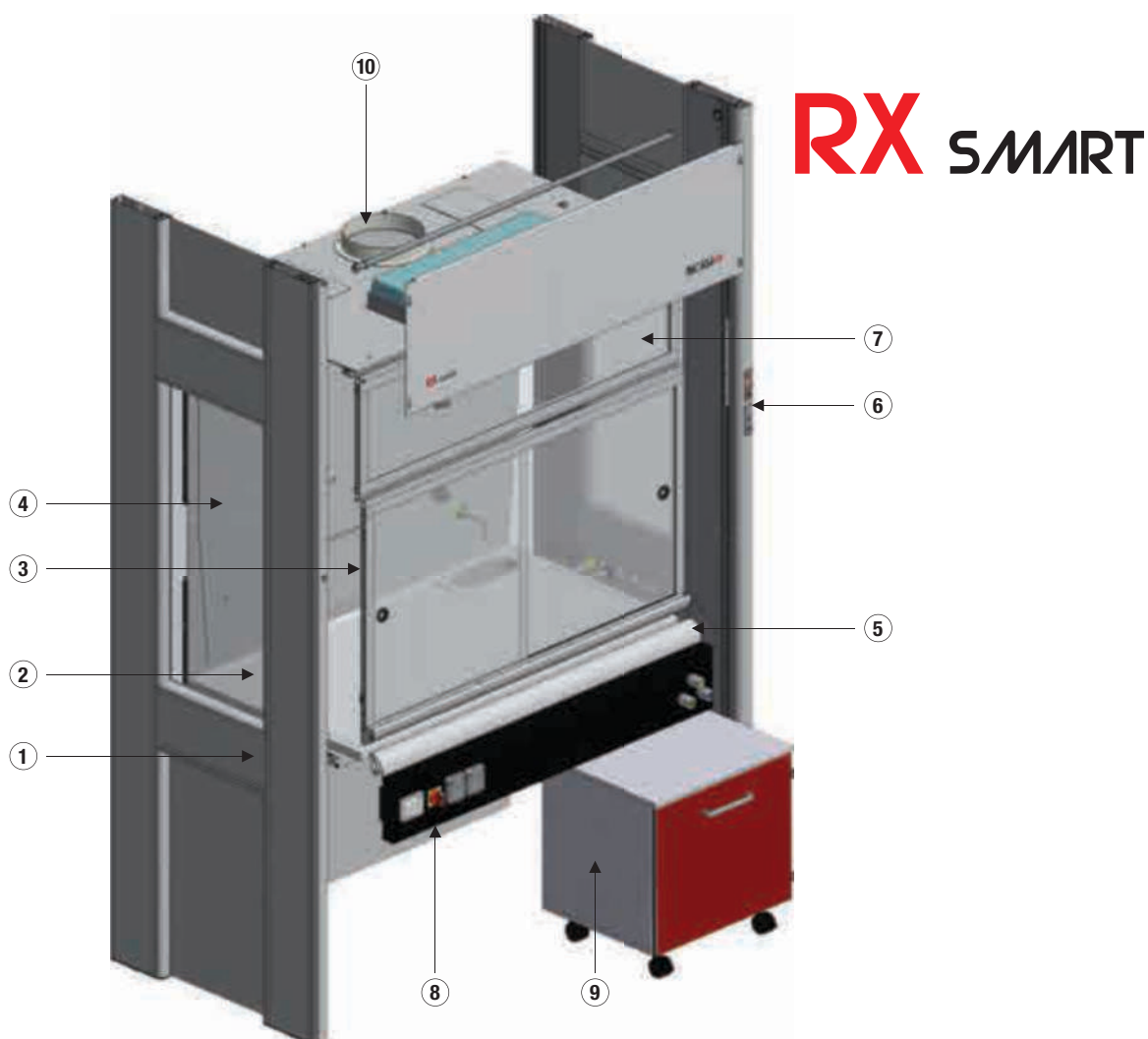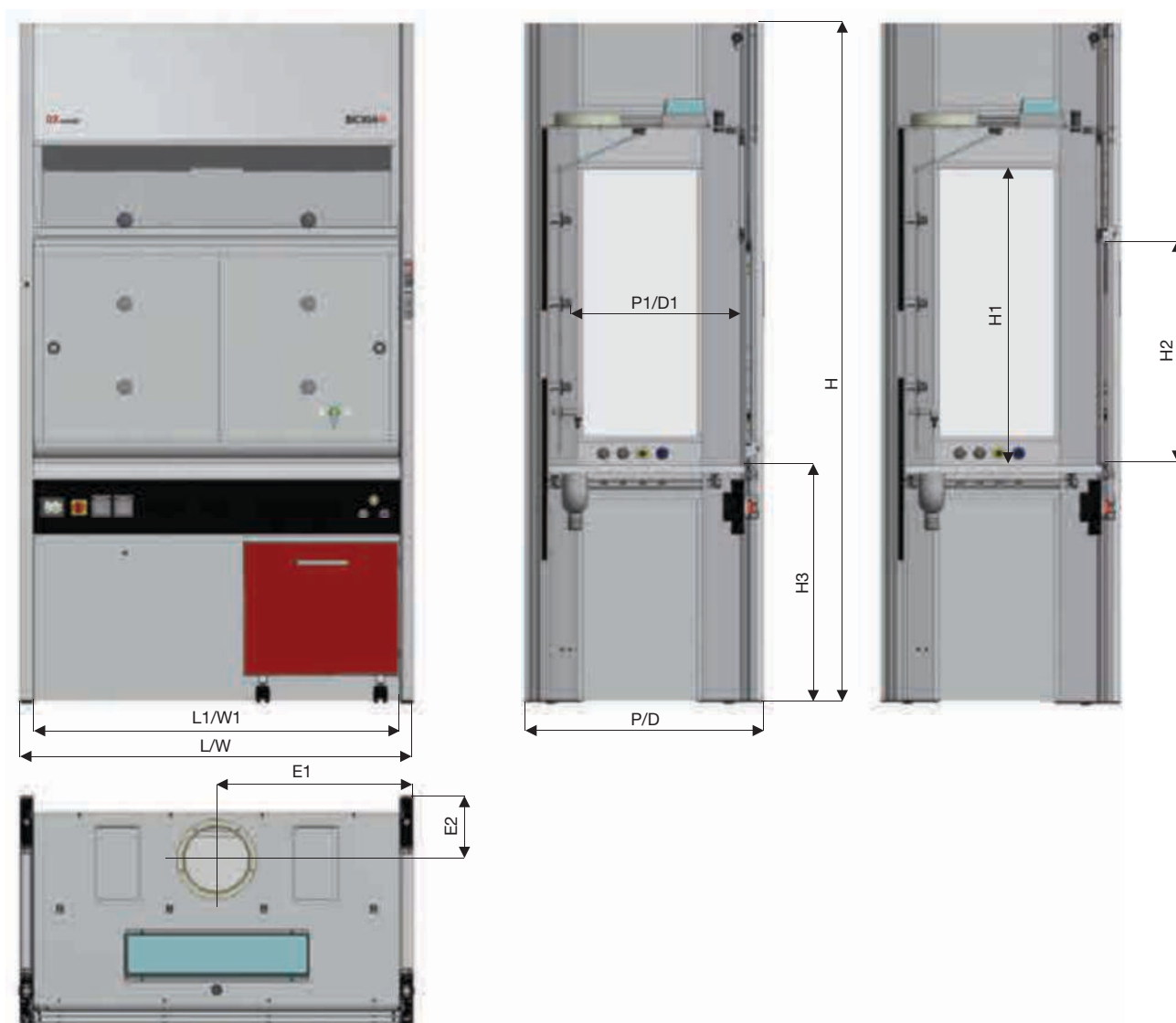
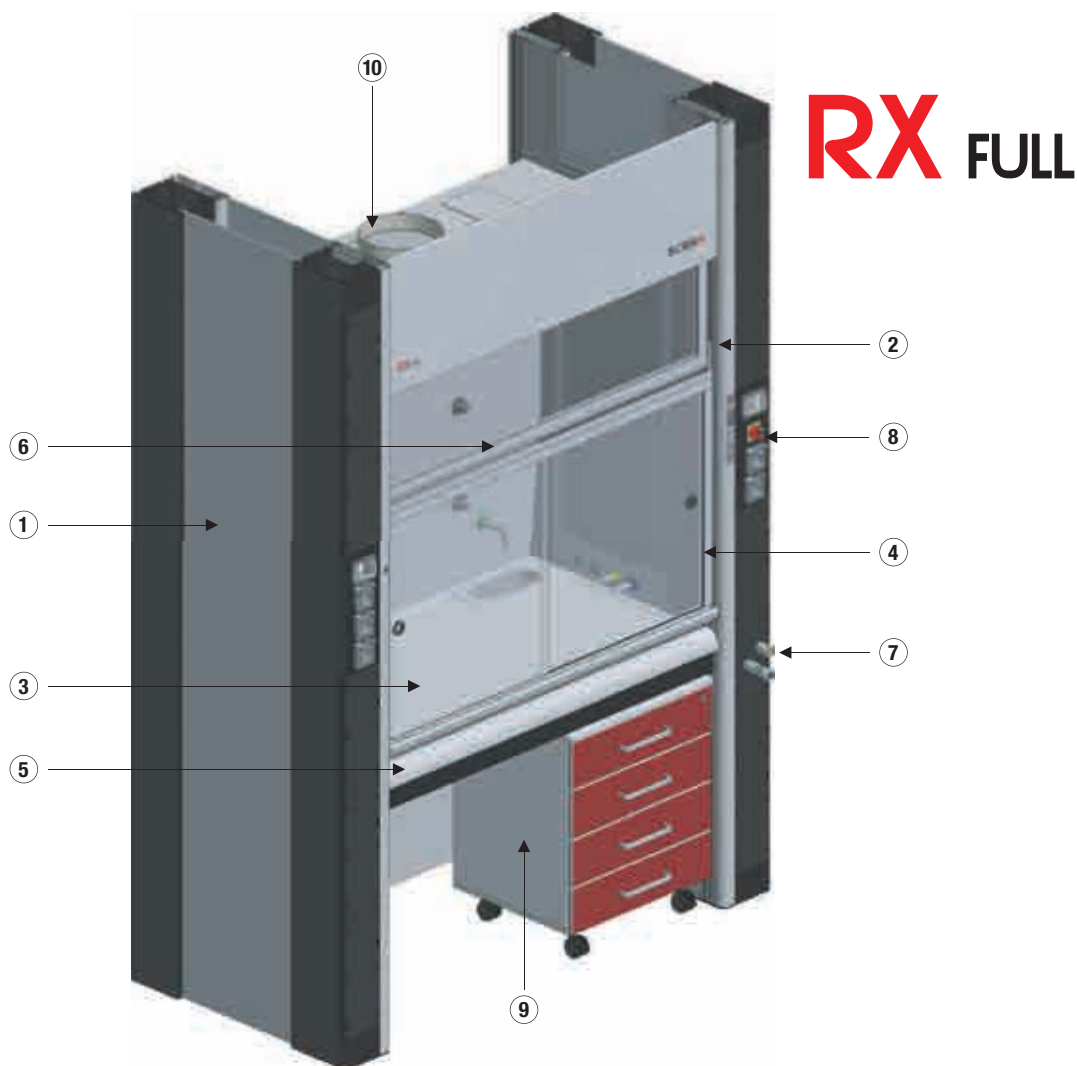
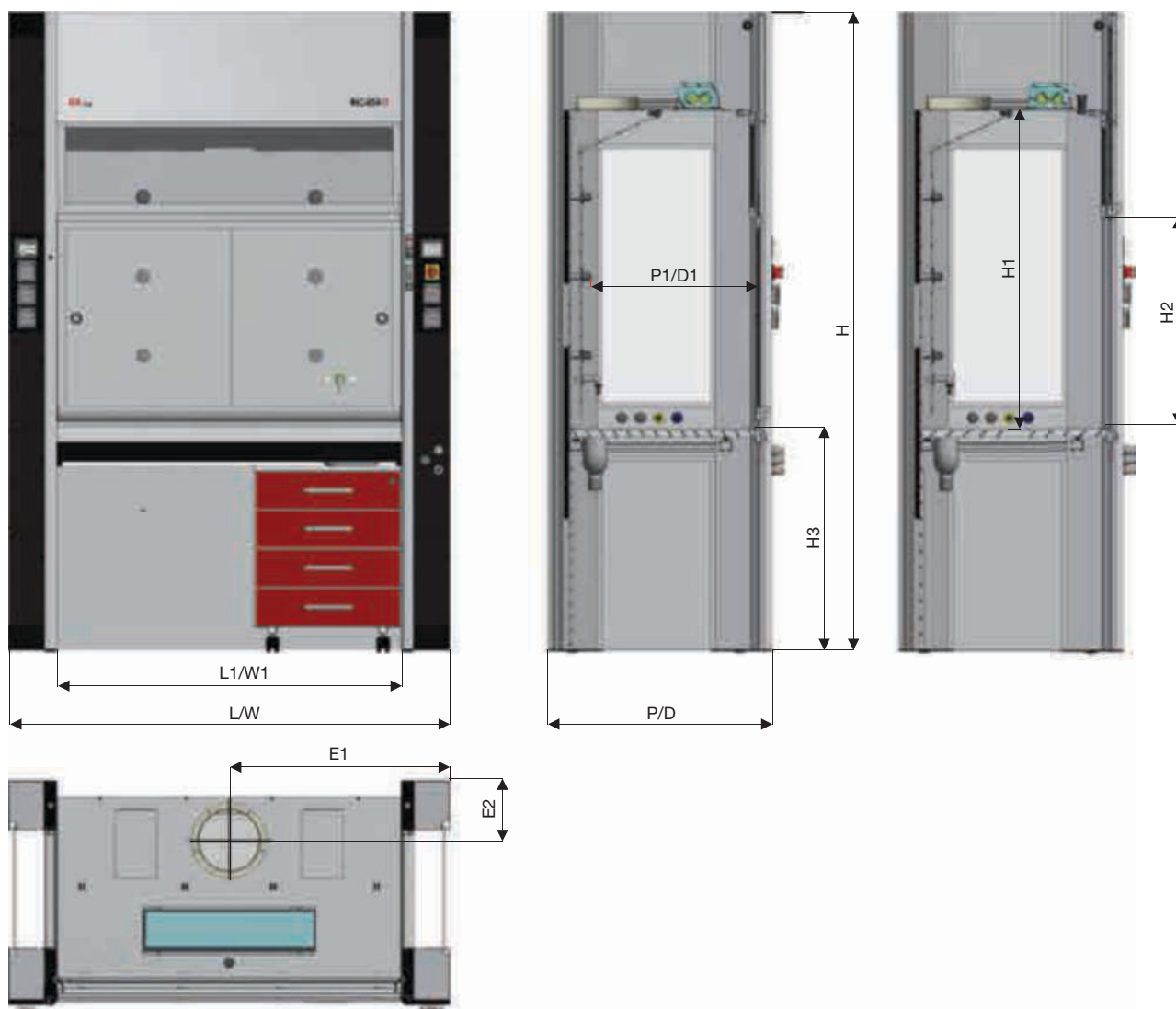
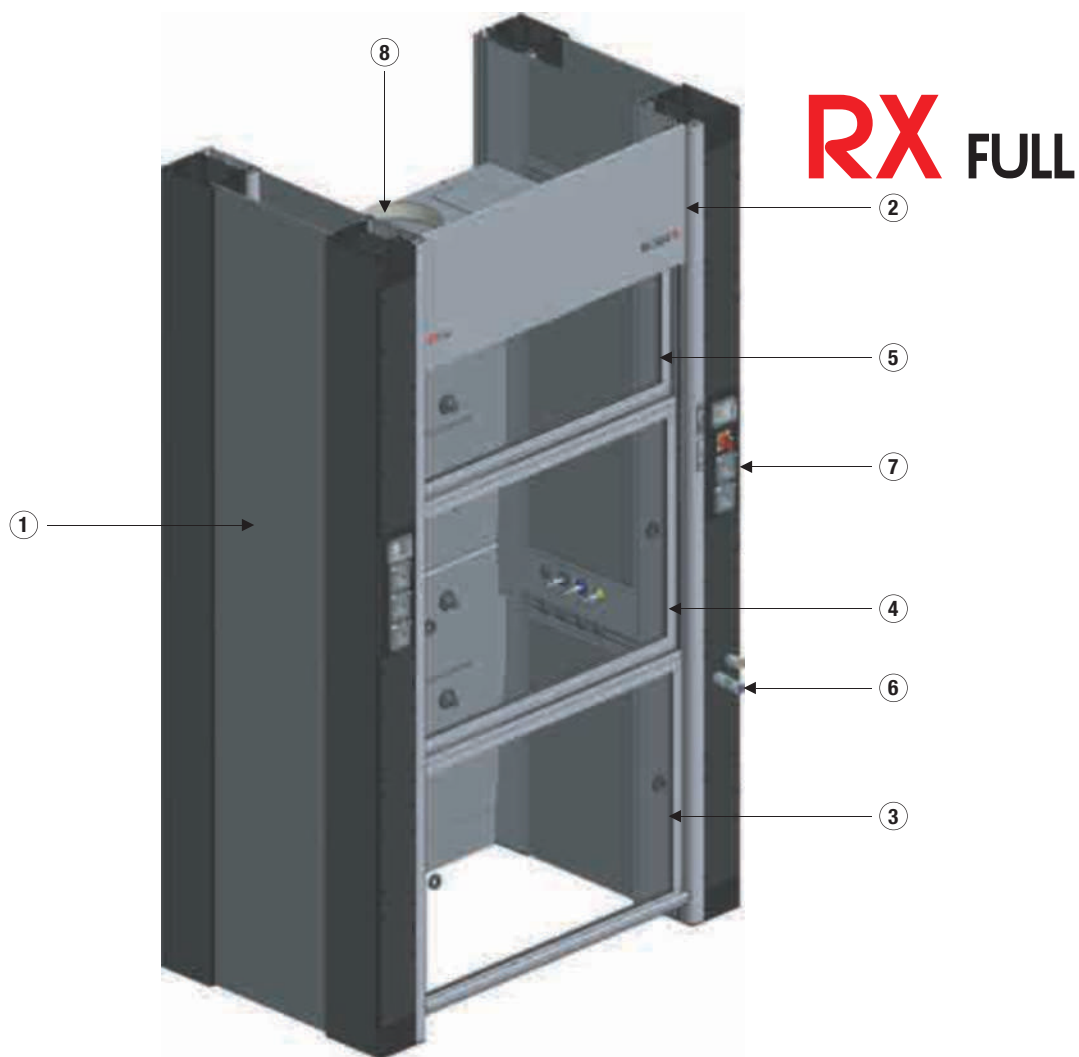
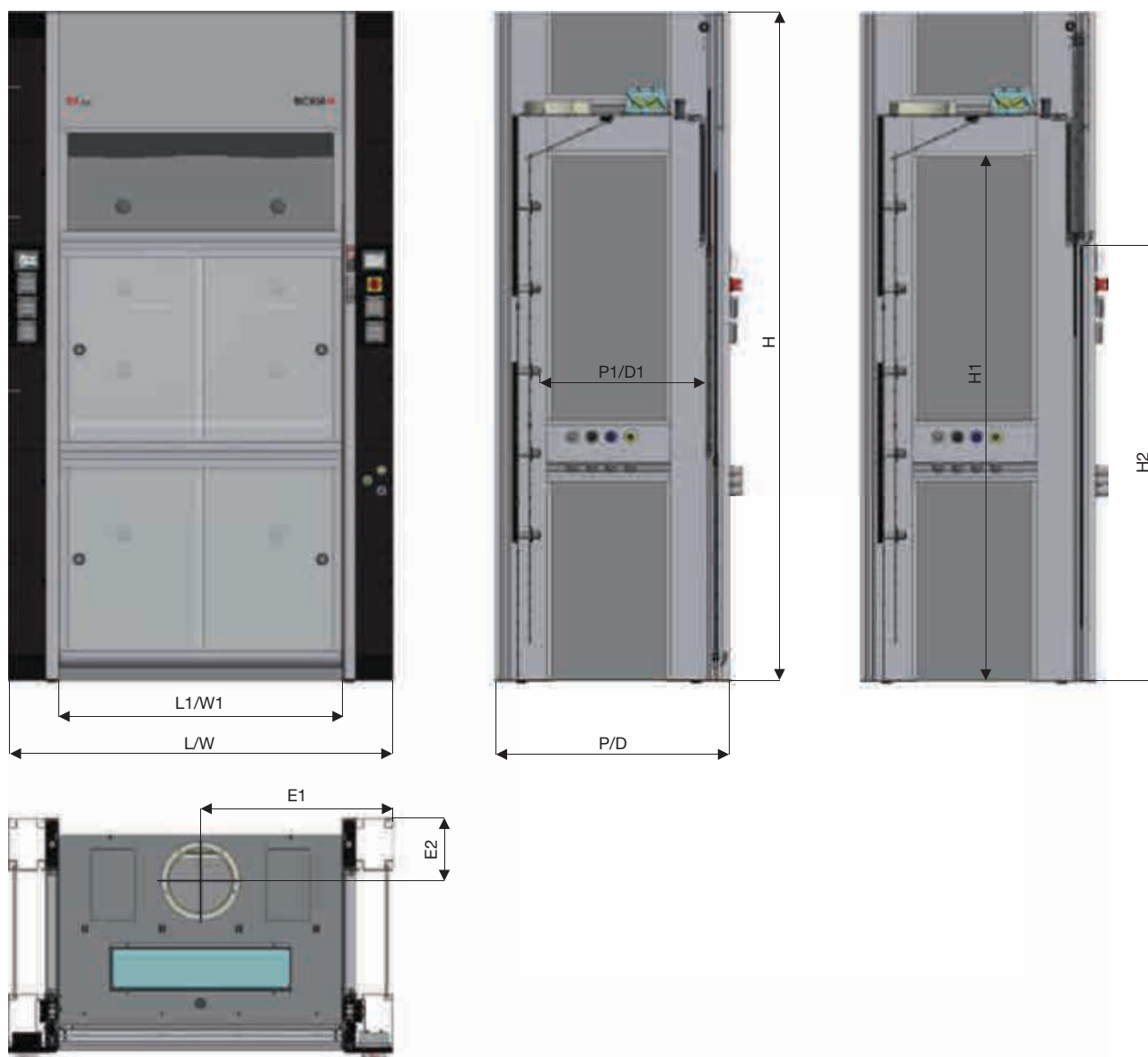
*A modern lab must provide maximum **integration among man, technology and environment**. Easy installation, flexibility, quality and safety are the essential guidelines leading Bicasa to the innovative design of IKARO, the functional supply system, and KOMPATTO, the dynamic furnishing system.*

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*Process automation, combined with advanced operative methods, draws new guidelines of laboratory design, suitable for workflow and instrumentation requirements. **IKARO suspended supply system** and **KOMPATTO mobile workstation** provide a perfect integration with the most advanced technologies with a complete flexibility in use.*

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*IKARO deletes forever the typical limits of traditional lab design. Built-in technologies applied to a suspended supply system simplify and optimize design, installation and testing phases of **modern lab**.*

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*IKARO supply beam gives a wide range of arrangements available to accommodate project and users requirements: starting from **FULL configuration**, the completely integrated solution, running to **SMART**, the slim type with small dimensions, ending to **WALL**, the wall mounted arrangement.*

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*IKARO has been designed as a modular system with the full integration of all components. The supply beam frame is composed of **modular units that can be combined to the lenght required**. Accessories and service outlet panels can be placed in the preferred position, easy to be changed or completed at any time.*

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*The rounded shape design is the distinctive characteristic of IKARO and KOMPATTO. The use of aluminium profiles increases the structural rigidity and reduces the overall weight of the suspended supply system. The snap fit joint system, used for aluminium profiles coupling, provides a direct access to the supply lines recesses.*

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
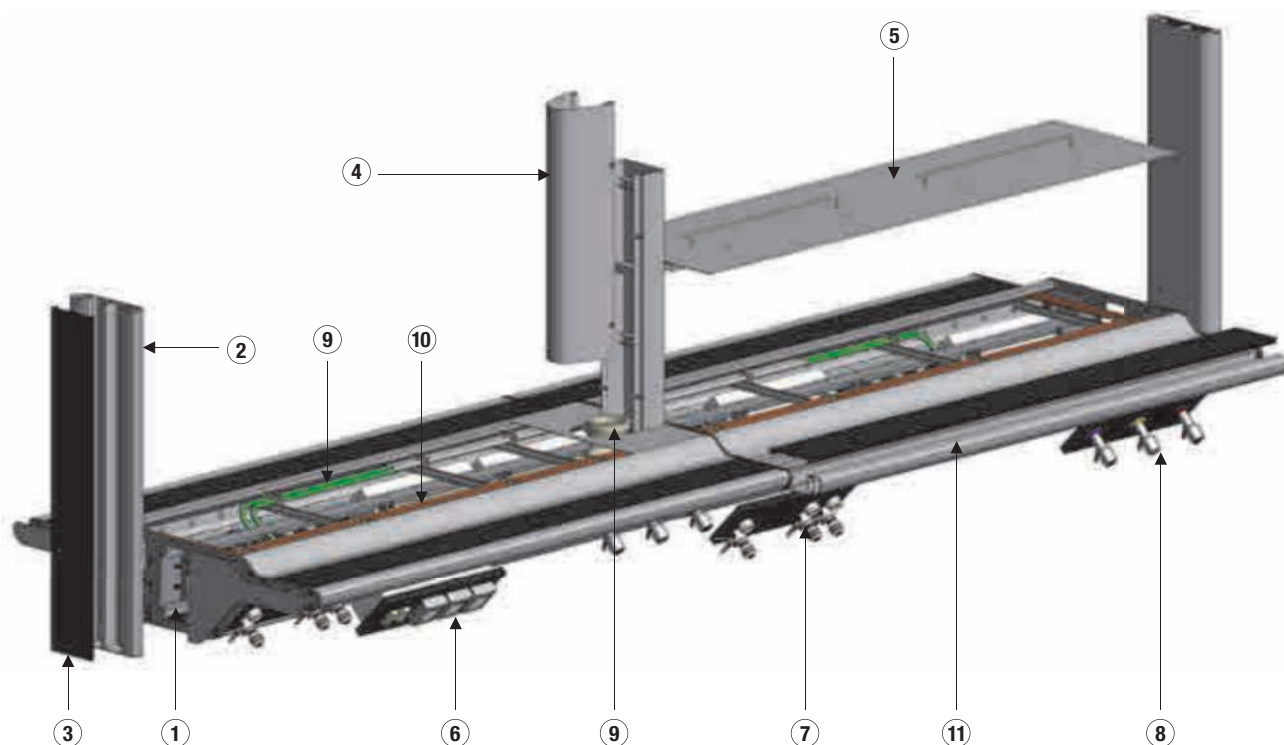
*IKARO supply beam has been **designed to be installed in any existing condition**. Vertical uprights can be anchored to ceiling or to floor using dedicated fixing plates that provide an easy and quick installation.*

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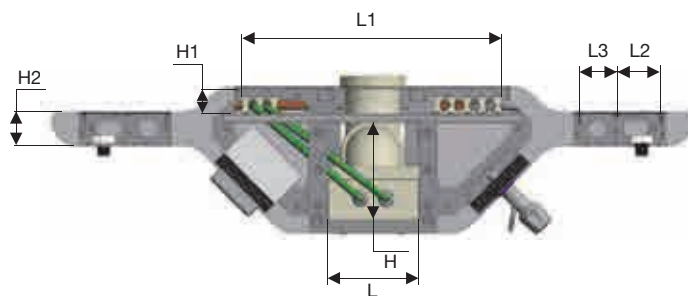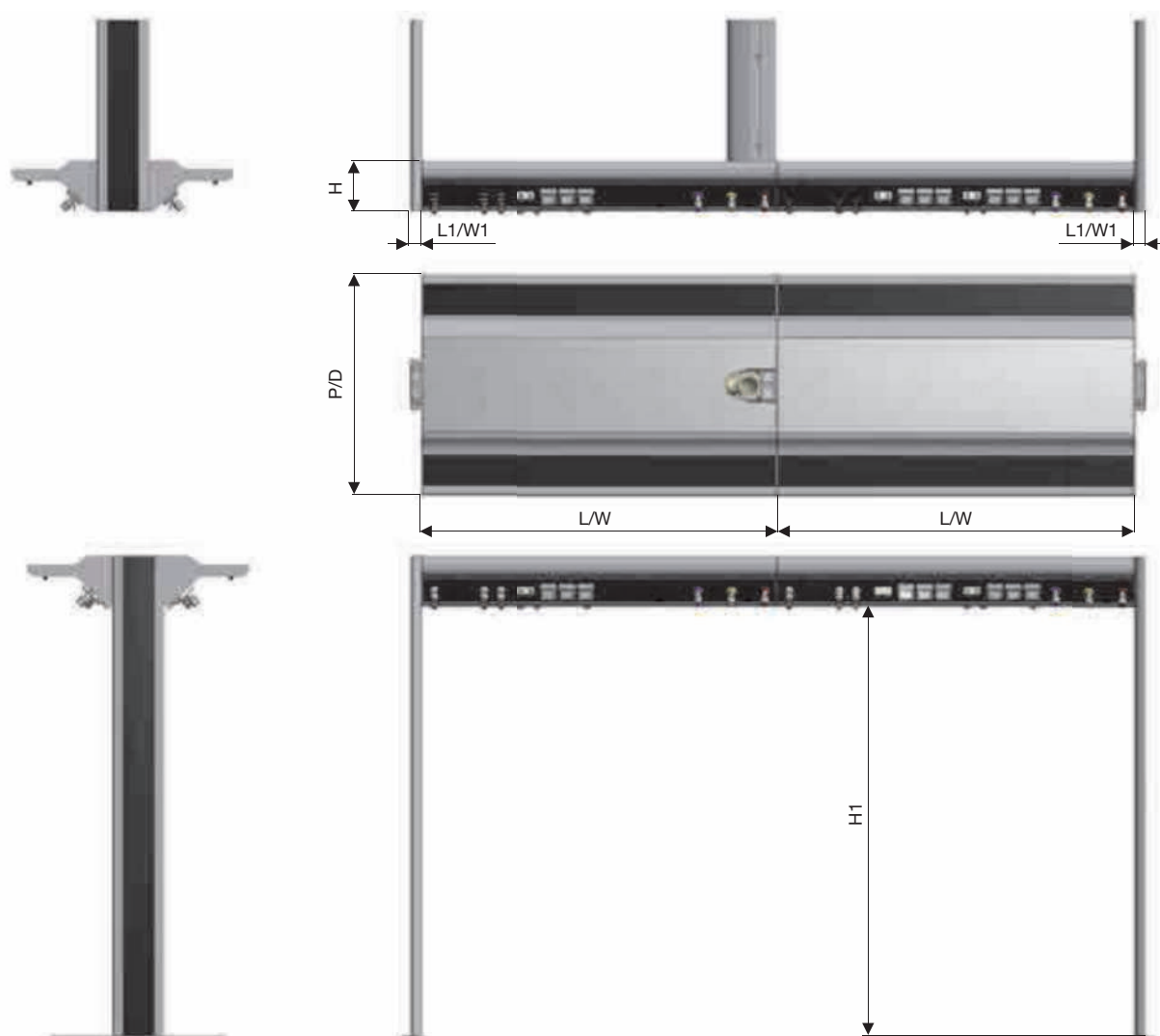
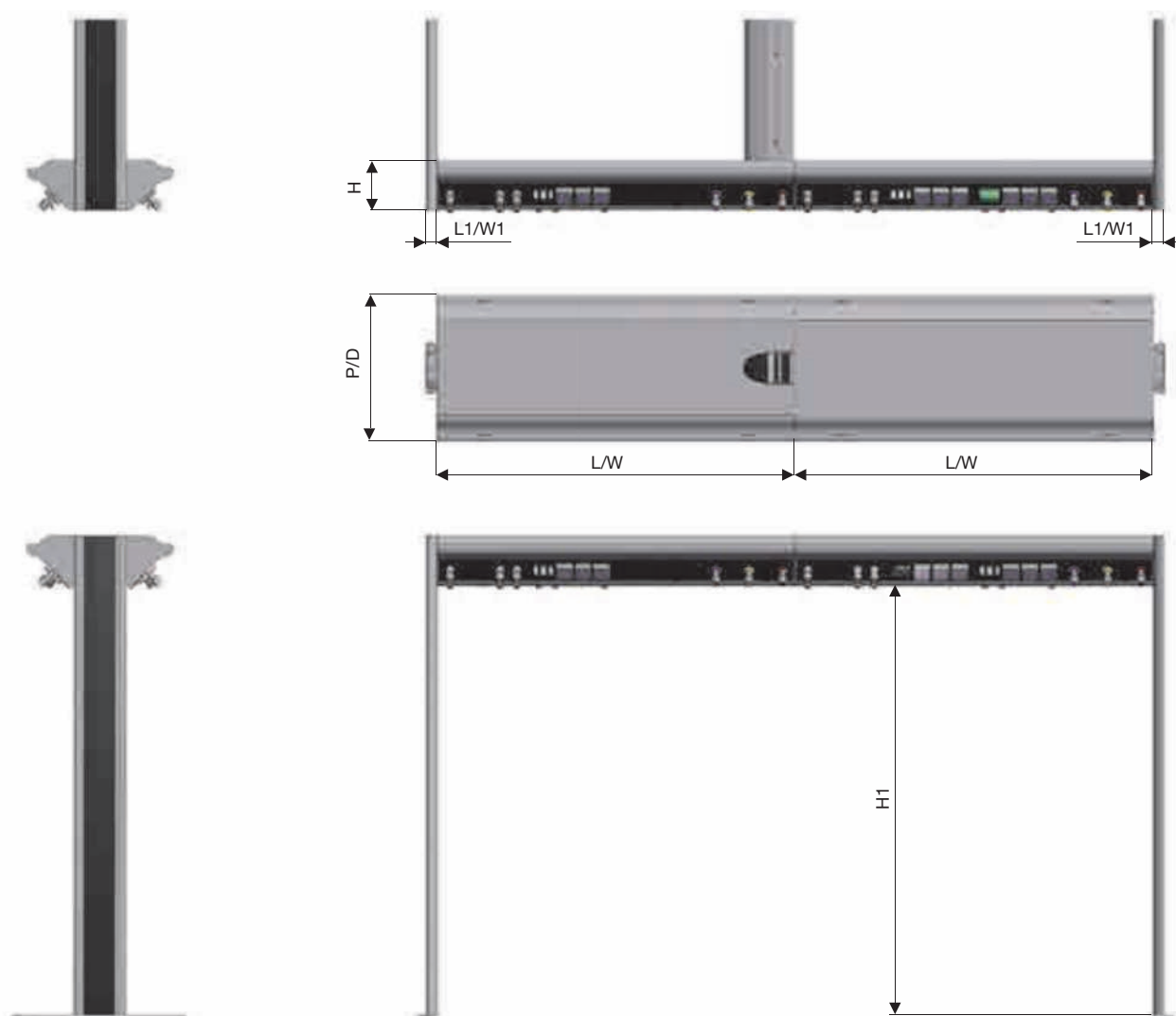
*The traditional lab changes face. The standard workstations change into KOMPATTO, **the dynamic high-tech system**. Steel and aluminum make KOMPATTO ergonomic, attractive and able to carry heavy loads.*

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*KOMPATTO mobile workstations are certified according to **EN 13150**, European standard concerning “Working benches for laboratories: dimensions, safety and test methods”.*

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*KOMPATTO mobile workstations are certified according to **BGR234/09.06** German standard, concerning workplace safety for storage devices. The load capacity tested and certified is up to 700 kgs, uniformly distributed, for all workstation types.*

## WORKSTATION PERSONALIZZABILI / CUSTOMIZABLE WORKSTATION

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*The dynamic and flexible system for ongoing needs. Starting from SMART, the basic configuration, running to FULL, the most complete solution, KOMPATTO fits with every laboratory. Shelves, monitor supports, CPU holder and cabinets increase the available space. Light bars and service outlets improve end user comfort.*

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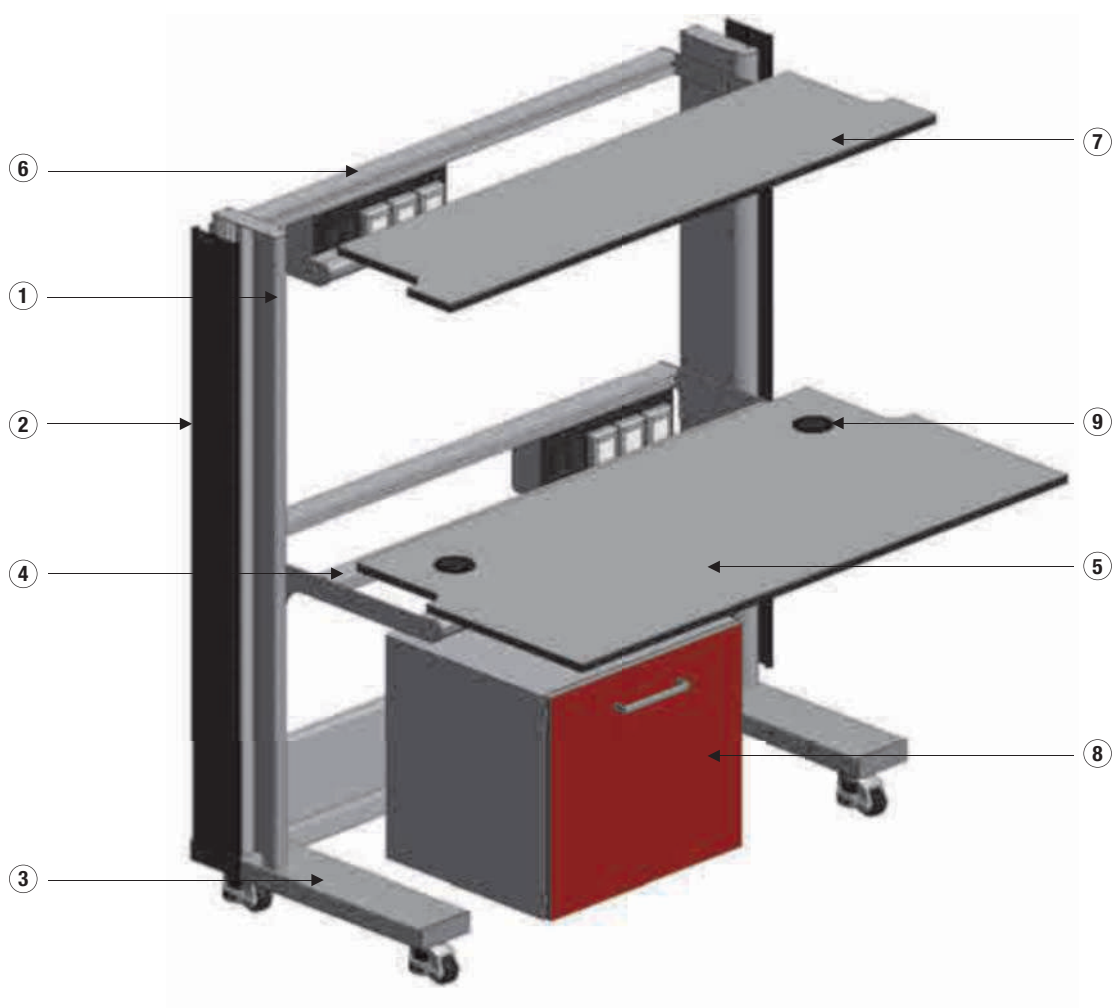
*KOMPATTO workstation lets you adapt the workplace to requirements, easily and quickly. Telescopic leveling castors, with brake system, change the mobile workstation into a fixed bench. Height adjustable countertop fits with every working requirement.*

## DATI TECNICI / TECHNICAL DATA

① Montante in alluminio  
*Aluminum upright*

② Carter rimovibile in alluminio  
*Aluminum removable panel*

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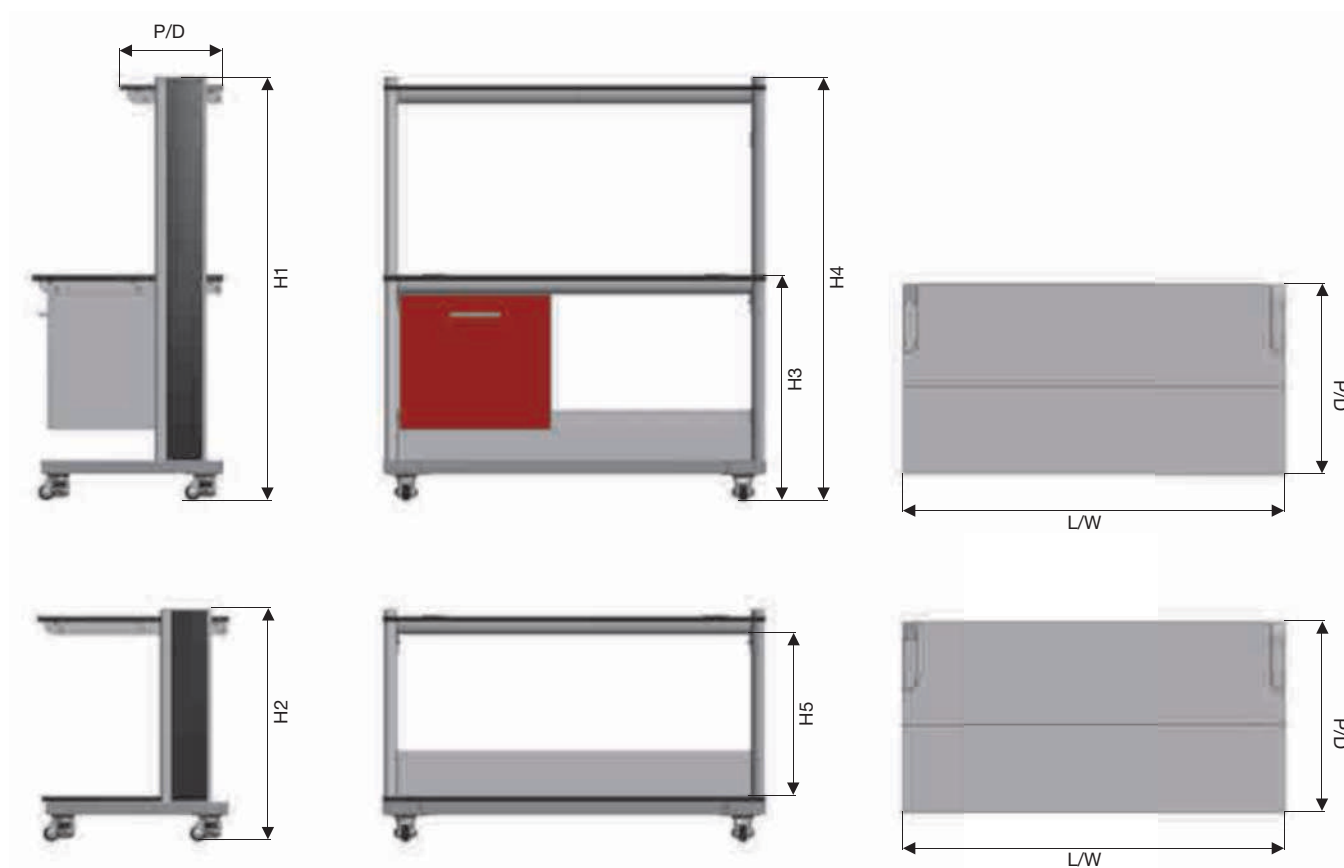
(\*) Il peso non include gli accessori e la mensola  
Weight value does not include accessories and overhead shelf.

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*COLUMNIA is the innovative laboratory supply system designed for functional and flexible point-of-use distribution with integrated piping and wiring. Suspended to media ceiling or to the concrete ceiling, the vertical column is best solution for projects requiring space as main task.*

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*COLUMNIA is designed to be factory assembled, testing the final configuration before the delivery with the pre-arranged pipes and wirings ready to be connected on site. Factory-tested process makes possible to decrease the duration of installation and commissioning phases with a relevant cost saving.*

## DESIGN E SICUREZZA / DESIGN AND SAFETY

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*The construction with a slim size and rounded shapes, including the HStrip panels ready to accept laboratory services, are the distinctive features of the vertical service column. Fast coupling system with snap-fit joints makes easier maintenance and reconfiguration after the installation. The access to the dedicated fittings compartment can be easily done both on the front and the back side of the column, removing single HStrip panels.*

## SOLUZIONI HI-TECH / HIGH-TECH SOLUTIONS

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*Framework is fully manufactured in aluminium - 100% recyclable material - consisting of Bicasa custom-made extruded profiles. The choice of such a material makes possible to decrease the weight drastically. The frame of the vertical column, designed to accept dedicated accessories, can be arranged with a huge quantity of service outlets, thanks to the HStrip horizontal panels easy to be customized as per the requirements.*

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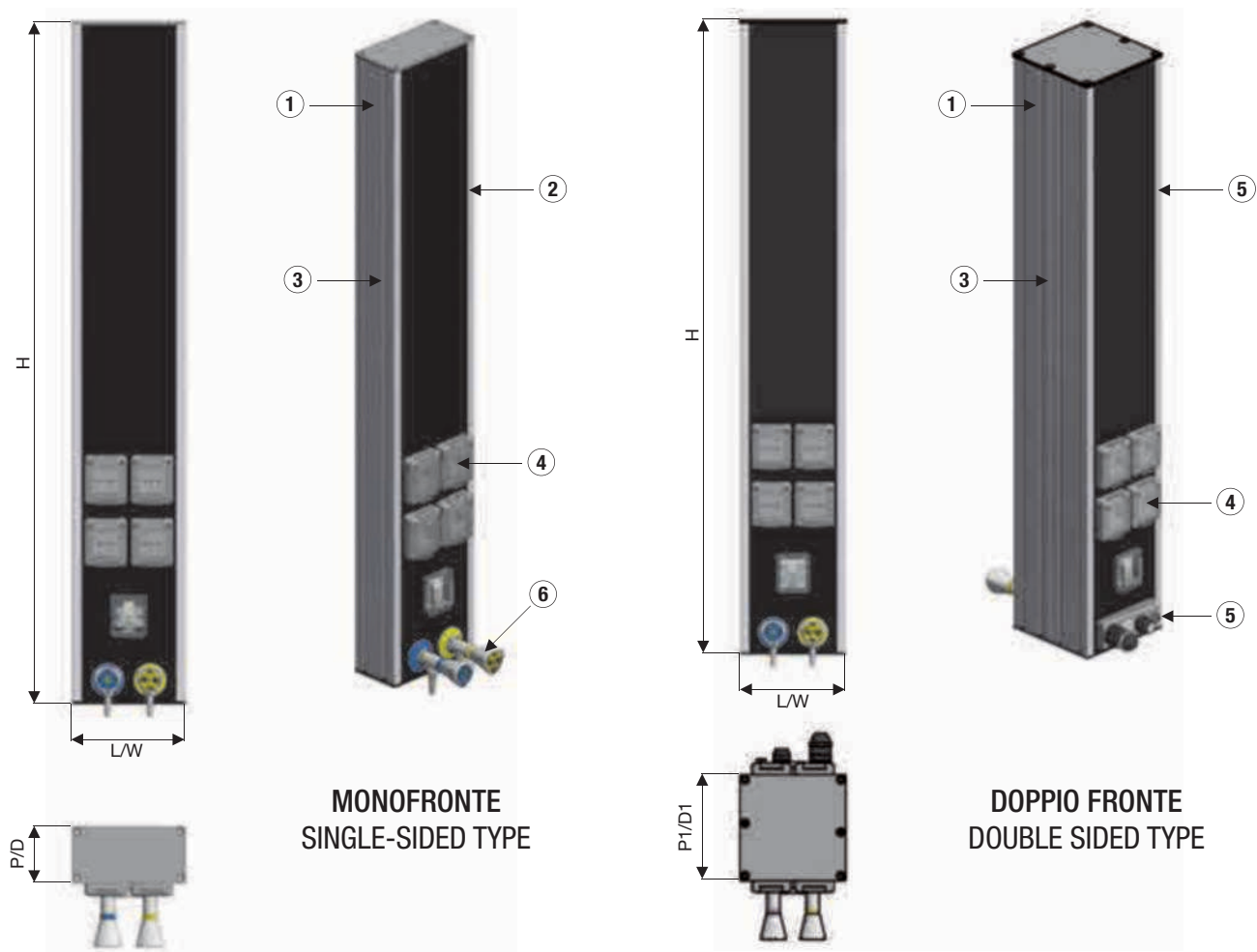
*The aluminium parts are treated with a polyurethane power coating suitable to protect surfaces and increase chemical resistance during laboratory use. The metallic matte light grey colour, consisting of Bicasa custom-made coating, is a remarkable innovation in aesthetic quality.*

# DATI TECNICI / TECHNICAL DATA

- ❶
 

Profilo in alluminio portante colonna  
*Vertical column main profile, aluminium made*
- ❷
 

Pannelli HStrip rimovibili porta servizi in alluminio  
*HStrip removable panels for service outlets, aluminium made*
- ❸
 

Binario guida per installazione di accessori sospesi (mensole, armadietti, ecc.)  
*T-slot track suitable for installation of suspended accessories (shelves, cabinets, etc.)*

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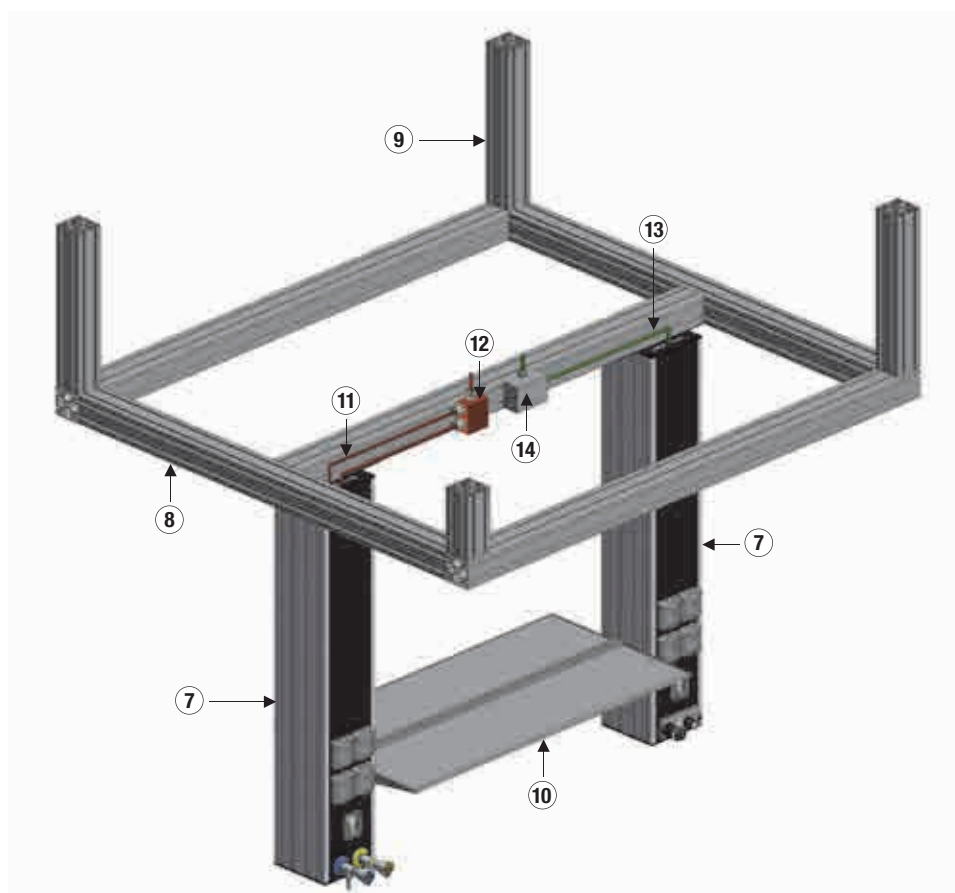
## GRIGLIA MEDIA CEILING / MEDIA CEILING CONCEPT

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*Laboratory projects require a perfect coordination between end user's requirements and the existing architectural features. Bicasa can **work in synergy with construction firms and architects** to achieve the highest functionality and the perfect integration, developing customized solutions for each project.*

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*KRISTAL is the exclusive furniture system designed, engineered and manufactured for projects in which **integration between laboratory casework and raised floor installation** is required. The self-supporting column is anchored below access flooring, ready to accept horizontal pipe lines and wiring.*

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*The vertical column, **fully aluminum made**, is designed to accept service outlets, minimizing the space required. Supply lines are arranged in vertical separated fittings raceways. Suspended countertops, under worktop cabinets and overhead accessories, are height adjustable.*

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*KRISTAL furniture system is designed ready to be integrated with partition elements, in order to arrange single working areas. The overall depth of 36 mm minimizes the space required for installation and can be equipped with different surfaces vertical panels.*

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*The flammable safety cabinets are designed for the storage of flammable materials in the working areas. They are certified in compliance with EN 14470-1 European standard, tested type 90 with a fire resistance of 90 minutes. Safety cabinets are ready to work with ventilation system for exhaust air.*

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*The chemical safety cabinets are designed for the storage of aggressive, non-flammable hazardous materials in working areas. The acids and alkalis cabinets consist of two hermetically separate storage chambers, with the possibility to connect a ventilation system for exhaust air.*

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*The construction of the safety storage cabinets increases the protection against corrosion and provides surfaces easier to be cleaned. Doors, including a cylinder locking system, can stop in any opening position. The base of safety storage cabinet is designed with integrated transport device for easy moving.*

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*Storage safety cabinets are available in a wide range of models with different sizes of width and height, depending on end user requirements. The internal configuration can be provided with high-adjustable shelves or with sliding drawers, heavy-duty type, for a safe and easy way of working in lab.*

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*Cupboards and underbench cabinets are designed for material storage. A wide range of choices allows to define **customized arrangements**, supplied with doors, drawers, glass panels and hinged or sliding opening. Elements can be added with horizontal and vertical integration, with optional rails and ladders to reach high placed shelves.*

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*Underbench cabinets and cupboard are certificated according to **EN-16121**, European standard concerning "Non-domestic storage furniture - Requirements for safety, strength, durability and stability".*

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*Cabinet carcass and top are manufactured in light grey colour. **Front panels of doors and drawers can be customized** as per end-user preferences. The wide range of colours available includes: maroon, charcoal grey, blue and light grey.*

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*Cabinet carcass and front are made of resin-bonded particleboard panels, fireproof type class B-s1, d0 according to EN 13501-1. **Bicasa works only with ecological panels, made of recycled wood.** Steel cabinet fabrication is available on request.*

*Each door or drawer is equipped with aluminum handle, easy to clean. Height-adjustable shelves and cabinet base are made of epoxy coated steel. Underbench cabinets are provided with swivel casters including front brake system.*

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*Drawers are assembled with steel sides height of 85 mm. The sliding system supplied ensures **self-closing** and load capacity up to 30 kgs. All drawer cabinets are provided with central lock and tilting-proof device. Extension sides, soft-close system and fast removing device of frontal panel are available on request.*

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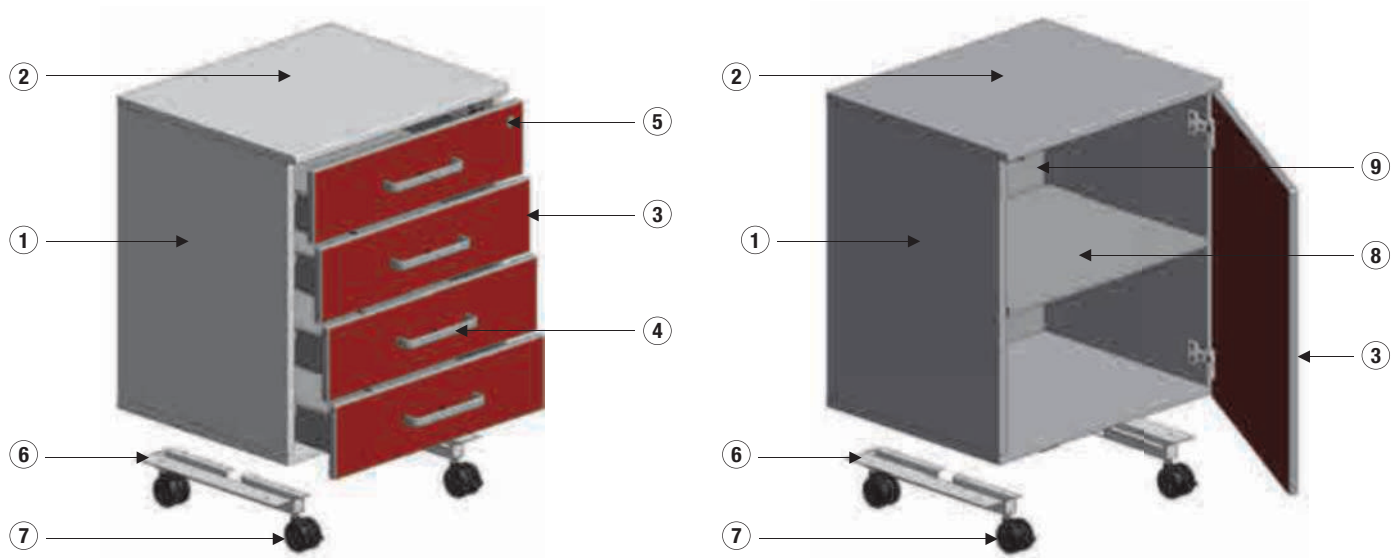
*Cabinets are **available with hinged or sliding doors**. Both the configurations can be equipped with melaminic panel or safety glass. Single pivot hinges work with 270° wide-angle opening. Concealed hinges are available on request. Closing profiles are supplied with dust-tight strip.*

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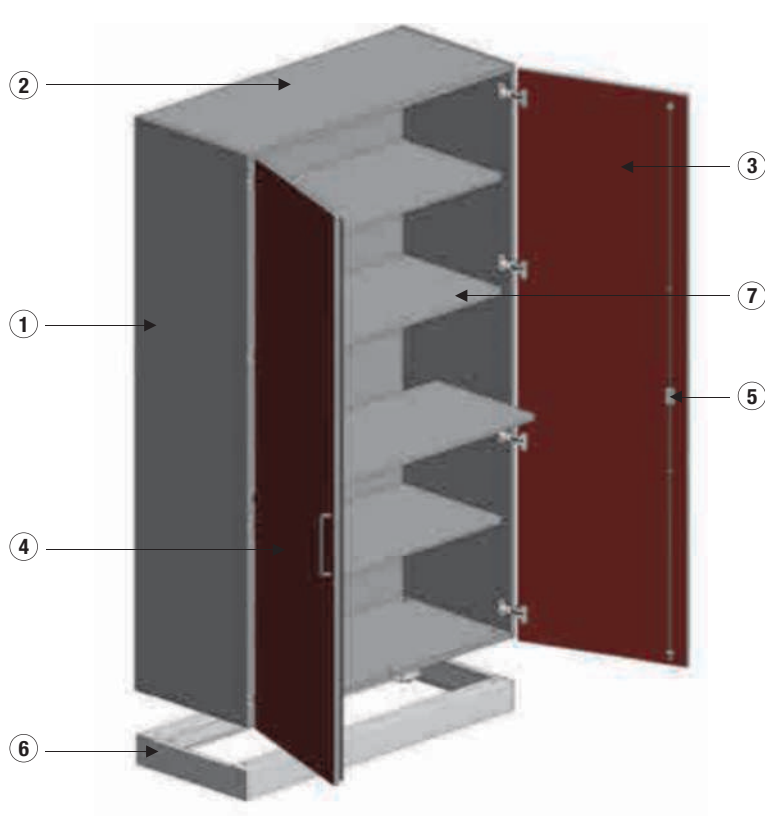
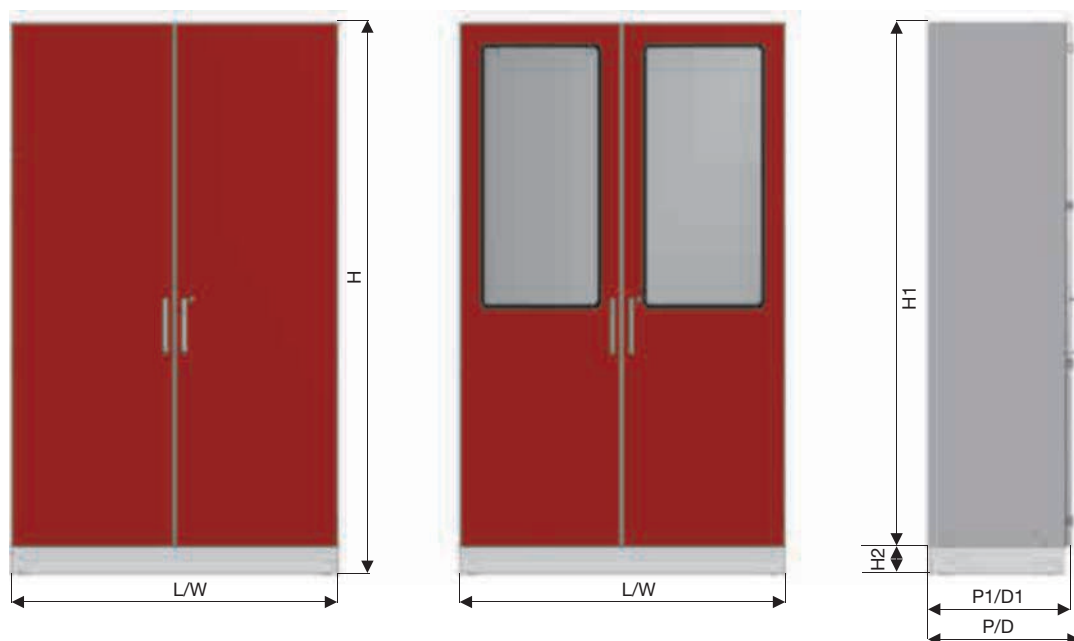
## MATERIALI / MATERIALS

### LAMINATO PLASTICO / PLASTIC LAMINATED

SPESSORE / THICKNESS mm 30

BORDO / BORDER Piatto / Flat

COLORE / COLOUR

GRIGIO CHIARO / LIGHT GREY

### LAMINATO PLASTICO POSTFORMING / POSTFORMING PLASTIC LAMINATED

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**BICASA™** all rights reserved
BICASA™ all rights reserved.